

## Use Case: Leveraging MagLev for Separating Powdered Illicit Drugs in Law Enforcement

### Introduction

Law enforcement agencies face challenges in quickly and accurately analyzing seized drug samples, which often contain complex mixtures of substances at varying concentrations. Traditional methods can be time-consuming, require sophisticated equipment, and may not be sensitive enough to detect low-concentration synthetic opioids like fentanyl. The MagLev device—using a novel approach called **magneto-Archimedes levitation**—offers a rapid, field-adaptable solution by separating drug mixtures based on density. This approach allows law enforcement to screen for multiple components in powdered samples, improve presumptive identification, and streamline lab analysis.

### How MagLev Technology Operates

MagLev technology functions by placing a cuvette filled with a paramagnetic solution (such as gadolinium(III) chelate in a non-polar solvent) between two magnets with their like poles facing each other. When a powdered sample is introduced:

- 1. Density-Based Separation:** The magnetic field causes compounds in the mixture to levitate at different heights based on their density. The precise levitation height correlates with each compound's density, allowing for visual stratification of multiple substances within the same sample.
- 2. Layered Analysis:** In practice, each levitating layer corresponds to a unique component or mixture of components in the drug sample. For instance, in a mixture of fentanyl, heroin, and lactose, each will float at a different level, creating distinct layers. This setup enables officers and lab technicians to separate even low-concentration active drugs from diluents and adulterants, which are typically added to increase volume or mimic specific properties.


- 3. Compatibility with Further Analysis:** After separation, individual fractions can be extracted with a pipette and subjected to confirmatory techniques like FTIR (Fourier-transform infrared spectroscopy) or Raman spectroscopy. This method improves the specificity and accuracy of subsequent analyses by isolating components beforehand.

### Key Applications in Law Enforcement Laboratories

**Enhanced Forensic Lab Analysis:** MagLev is designed to streamline forensic workflows by isolating compounds within complex drug mixtures based on density. In-house lab technicians can separate substances for more specific analysis, improving the accuracy of subsequent confirmatory methods such as FTIR and Raman spectroscopy. This is especially valuable for detecting low-concentration substances, like fentanyl, that are difficult to analyze directly in mixtures.

Learn More:  
[www.verichemtech.com](http://www.verichemtech.com)



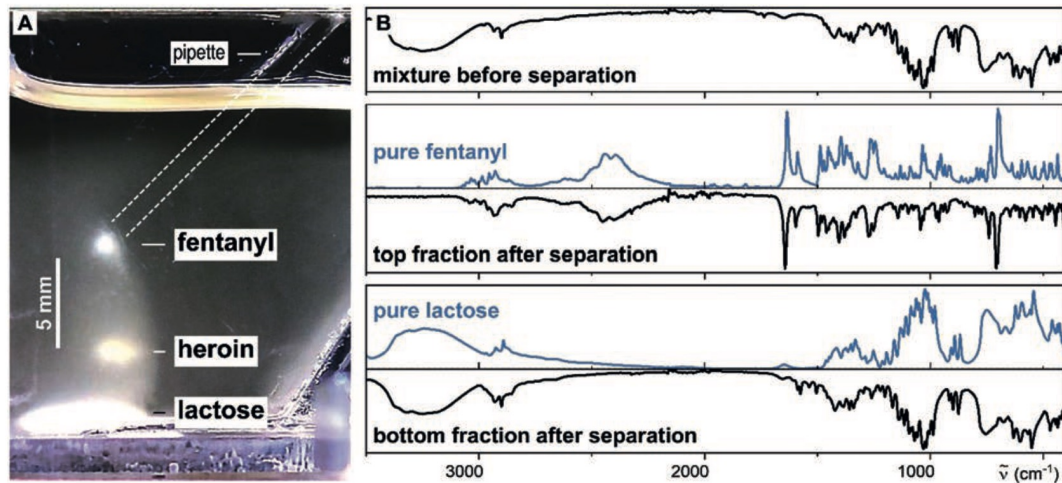

**Figure 1.** Successful separation, presumptive identification, and confirmatory identification of dilute fentanyl, in a mixture of heroin and a diluent (lactose). A) An image taken after 30 min of separation by MagLev of a powdered mixture of fentanyl-containing heroin (fentanyl-HCl (1.3 wt.%), heroin-HCl (2.6 wt.%), and a-lactose (96.1 wt.%)). The separation was performed in a custom-made cuvette (shaped to allow easy entry of a pipette) filled with a paramagnetic solution of Gd(DPM)3TOPO (450 mM) in a mixture of 23 vol% hexane and 77 vol% tetrachloroethylene. The separated fractions were extracted using a Pasteur pipette and were subsequently rinsed with hexane under suction filtration to remove any remaining gadolinium complex and air-dried. FTIR-ATR spectra (normalized to the highest peak) were measured from the powdered mixture before separation (top spectrum). The extracted fractions containing fentanyl and lactose (lower two black lines), and the pure compounds (blue lines) are below.

**Backlog Reduction and Faster Turnaround:** Forensic labs often face backlogs when dealing with intricate drug samples that require detailed analysis. MagLev provides a quick, efficient method for initial compound separation, enabling labs to process samples faster and more accurately. By isolating components upfront, MagLev allows for prioritization and quicker throughput, ultimately supporting a higher volume of sample processing.

**Improved Resource Efficiency:** With MagLev, forensic labs can conduct preliminary separations that reduce reliance on more labor-intensive methods. This preparatory step not only conserves resources but also enhances lab productivity by enabling faster, more targeted follow-up testing.

## Practical Benefits of MagLev Technology in Forensic Laboratories

MagLev provides essential advantages for forensic labs managing complex drug cases:

**Non-Destructive Testing:** Unlike traditional methods that consume the sample, MagLev preserves the material, enabling further confirmatory testing without compromising evidence integrity. This non-destructive approach ensures that each component in a sample remains available for thorough analysis.

**Reduced False Positives and Negatives:** Complex drug mixtures often lead to inaccurate results due to masking agents or interfering substances. MagLev reduces this risk by isolating individual compounds, minimizing the chance of false positives (e.g., legal substances flagged as drugs) or false negatives (e.g., potent drugs undetected due to dilution).

**Enhanced Laboratory Safety:** Forensic labs frequently encounter hazardous substances like fentanyl. MagLev's density-based separation occurs within a controlled solution, lowering the risk of accidental exposure and improving laboratory safety protocols for handling potent compounds.

**User-Friendly Operation:** MagLev's straightforward design requires minimal training, making it easy for lab technicians to incorporate into their workflow. The simplicity of its setup and clear visual separation of compounds enhances lab efficiency without the need for specialized technical skills.

**High Sensitivity for Diluted Samples:** Synthetic opioids, like fentanyl, are often highly diluted, complicating detection in mixtures. MagLev's ability to levitate and isolate even low-concentration substances (below 5%) enables reliable identification of these challenging compounds, providing forensic labs with a sensitive and accurate tool for complex cases.

## Conclusion

MagLev represents a significant advancement in laboratory-based drug analysis, offering an efficient, accurate, and safe method for handling complex drug mixtures. By isolating substances through density-based separation, MagLev enhances the effectiveness of confirmatory methods like Raman and FTIR spectroscopy, supporting forensic labs with a versatile tool to address the challenges of today's drug landscape. This technology improves workflow efficiency, increases sample throughput, and delivers high-quality results critical for modern forensic analysis.

Source: Adapted from "Analysis of Powders Containing Illicit Drugs Using Magnetic Levitation" by Christoffer K. Abrahamsson et al., *Angewandte Chemie International Edition* (2020). DOI: 10.1002/anie.201910177